

RNAZOL® BD PROTOCOL FOR THE EXTRACTION OF DNA FROM PLASMA OR SERUM.

William Wilfinger, Molecular Research Center, Inc.

The search for physiologically significant markers has spurred interest in the presence of extracellular DNA in circulating plasma or serum. Although RNAzol® BD is designed for the extraction of RNA from blood, this protocol employs alkaline reagent that allows isolation of minute quantities of DNA from plasma or serum samples.

LYSIS

1. Add 1 ml of either plasma or serum to 2 ml of RNAzol® BD.
2. Make sure the tube is tightly closed and vigorously shake the sample - RNAzol® BD mixture for 30 seconds.
2. The lysate may be processed immediately or stored at -20 C prior to DNA extraction.

PHASE SEPARATION

1. Add 50 µl of 6 M NaOH for each ml of plasma or serum in the lysate and mix.
If samples were frozen, remove the lysate from the freezer and add 50 µl of 6 M NaOH for each ml of plasma or serum in the lysate prior to thawing. Place the tubes in a warm water bath and thaw the contents with shaking until completely thawed.
2. Store the solution at room temperature for 5 minutes.
3. Add 300 µl of BCP (1-bromo-3-chloropropane, MRC Catalog # BP 151) for each ml of plasma or serum in the lysate. Shake vigorously for 30 seconds and store at room temperature for 15 minutes.
4. Centrifuge at 12,000 g for 15 minutes at room temperature.

DNA PRECIPITATION

1. Transfer the clear water supernatant to a clean tube, add 1 - 2 µl of Polyacryl Carrier (MRC Catalog # PC 152) and mix.
2. Add an equal volume of isopropanol and thoroughly mix the solution. Store the sample for 30 minutes at 4 C.
3. Centrifuge at 12,000 g for 15 minutes at room temperature to obtain a pellet with DNA.

DNA WASH

1. Decant or remove the supernatant from the tube without dislodging the DNA pellet. Remove residual solution with a clean pipet tip.
2. Add 1 ml of 75% ethanol and vortex the solution. Store for 2 - 3 minutes at room temperature and centrifuge at 12,000 g for 1 - 2 minutes at room temperature. Decant or remove the wash solution with a pipet tip.
3. Perform a second wash with a fresh aliquot of 75% ethanol and remove all residual wash solution.
4. Centrifuge the pellet for 20 seconds at 12,000 g and carefully remove any residual wash solution.

DNA SOLUBILIZATION

1. Add 20 µl of water and solubilize the DNA pellet for 5 - 10 minutes with periodic vortexing.
2. The expected DNA recovery is below the detection limit of the Nanodrop spectrophotometer. Typically, 2 - 4 ul of the undiluted DNA solution used for PCR yields Ct values ranging from 20 - 40 depending on DNA copy number and primer efficiency and specificity.